

# MCCP Program (Multiple Column Cleaning & Pre-Conditioning) Copyright 2013

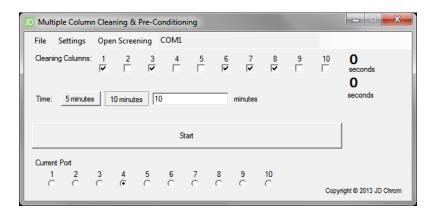

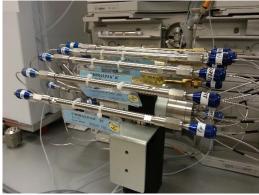

## ITEM DESCRIPTION

Included in the MCCPS System, the MCCP Program is integral to Phase 1 (Cleaning & Pre-Conditioning) of the MCCPS System's specialized 2-Phase Method Development Process. The streamlined user interface provides a fast, easy, and intuitive way to set up a cleaning & pre-conditioning process for upwards of 10 columns.

## **KEY FEATURES**

- Automated Column Cycling/Selection
- Resource Savings via Specialization
  - o Time, Samples, Solvents
- Simple & Intuitive User Interface
- Easy Column Selection via MCCP Program UI
- Time Display
  - Current Column & Total Run Time

## MCCP PROGRAM'S ROLE IN SPECIALIZED 2-PHASE METHOD DEVELOPMENT

To illustrate, let us assume the following in this example: 10 unique columns; 5min/column Cleaning & Pre-Conditioning Method; 20min/column Screening Method (injection); Let D = Distraction Cost (for user to put down everything & return to the HPLC System). Also note a common practice of injecting a sample 3x to achieve: (1) column equilibration (2) sample injection (3) injection confirmation.

### **Conventional Method Development**

Time	Task		
2	Column Setup		
63	Sample Analysis		
	1	Configure Method	
	2	System Startup (reaching method parameters)	
	20	Injection - C&P	
	20	Injection - Analysis	
	20	Injection - Analysis Confirmation	
1	Column Removal		
D	Distraction Cost		
66+D	Column Total		
	x(10)		
	FINAL TOTAL		
	660min+10D = 11hrs + 10(Distraction Cost)		

MCCPS System Method Development

Time	Task		
20	Column Setup		
		2min x 10columns	
53+D	Column Cleaning & Pre-Conditioning		
	1	Configure C&P Run	
	2	System Startup	
	50	Specialized C&P (5min x 10columns)	
	D	Distraction Cost	
203+D	Column Screening		
	1	Configure Screening Run	
	2	System Startup	
	200	Specialized Screening (20min x 10columns)	
	D	Distraction Cost	
276+2D	Total		
	FINALTOTAL		
	276min+2D = 4hrs 36min + 2(Distraction Cost)		

Total Time Saved in PHASE-1: 3 hours and 27 minutes & 10 times fewer Distraction Costs

260 minutes & 10 Distraction Costs VS 53 minutes & 1 Distraction Cost



PO Box 19146 Email: <a href="mailto:support@jdchrom.com">support@jdchrom.com</a>
Newbury Park, CA 91319-9146 Website: <a href="mailto:www.jdchrom.com">www.jdchrom.com</a>